

## DNA purification products from MACHERY-NAGEL

**BIOKÉ**  
sharing knowledge

Schuttersveld 2  
2316 ZA Leiden  
The Netherlands  
T. +31 (0)71 7200 220  
T. Belgium: 0800 71640  
F. +31 (0)71 8910 019  
info@bioke.com  
www.bioke.com

*DNA mini kit for microorganisms*

*Get microbial DNA the easy way*

**NucleoSpin<sup>®</sup> Microbial DNA**



**Suitable for a large variety of starting materials**

High quality DNA from gram-positive and gram-negative bacteria, yeast, and fungi

**NucleoSpin<sup>®</sup> Bead Tubes for efficient lysis included**

Compatible with many disruption devices

**Fast and easy procedure**

Liquid Proteinase K included – no additional enzyme required

**...high quality DNA for reliable downstream applications**

**MACHERY-NAGEL**

[www.mn-net.com](http://www.mn-net.com)




# NucleoSpin® Microbial DNA


## Product at a glance

Technology	Silica-membrane technology
Sample material	< 40 mg wet weight cell pellet
Fragment size	200 bp–approx. 50 kbp
Typical recovery	Depends on sample type and disruption Approx. 5–25 µg (30 mg wet weight cell pellet)
A <sub>260</sub> /A <sub>280</sub>	1.6–2.0
Elution volume	100–200 µL
Preparation time	35 min/6 preps
Binding capacity	60 µg

## Procedure

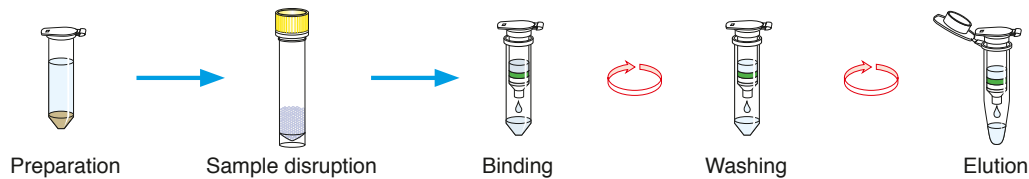

## Application data

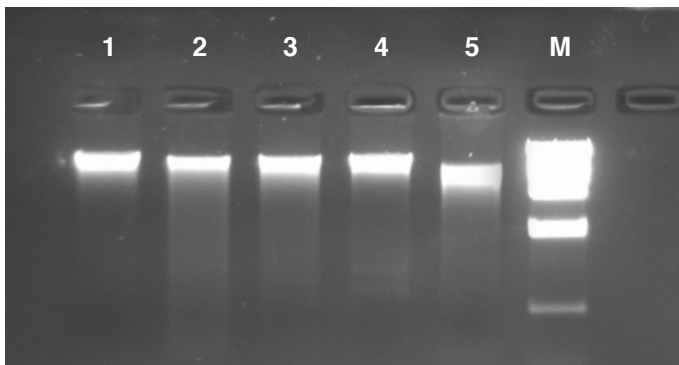

### Efficient DNA recovery from different microorganisms

DNA was isolated with the NucleoSpin® Microbial DNA kit and NucleoSpin® Bead Tube Type B (included in the kit) or NucleoSpin® Bead Tube Type C (see ordering information). 100 ng DNA per prep was analyzed by agarose gel electrophoresis showing high molecular DNA without RNA contamination or DNA degradation.

1. *Escherichia coli*, NucleoSpin® Bead Tube Type B
2. *Vibrio fischerii*, NucleoSpin® Bead Tube Type B
3. *Bacillus subtilis*, NucleoSpin® Bead Tube Type B
4. *Corynebacterium glutamicum*, NucleoSpin® Bead Tube Type B
5. *Saccharomyces cerevisiae*, NucleoSpin® Bead Tube Type C

## Ordering information

Product	Specifications	Preps	REF
NucleoSpin® Microbial DNA	NucleoSpin® Bead Tubes Type B included	10/50	740235.10/.50
Related products	Specifications	Preps	REF
NucleoSpin® Bead Tubes Type A	0.6–0.8 mm ceramic beads	50	740786.50
NucleoSpin® Bead Tubes Type B	40–400 µm glass beads	50	740812.50
NucleoSpin® Bead Tubes Type C	1–3 mm corundum	50	740813.50

### Microorganism

Microorganism	Tested by
<i>Acinetobacter</i>	customer
<i>Aspergillus spec.</i>	MN
<i>Bacillus subtilis</i>	MN
<i>Clostridium ljungdahlii</i>	customer
<i>Corynebacterium glutamicum</i>	MN
<i>Escherichia coli</i>	MN
<i>Eurotium spec.</i>	customer
<i>Klebsiella pneumoniae</i>	customer
<i>Microbacterium spec.</i>	customer
<i>Pichia pastoris</i>	customer
<i>Pseudomonas aeruginosa</i>	customer
<i>Rhizopus spec.</i>	MN
<i>Saccharomyces cerevisiae</i>	MN
<i>Staphylococcus epidermis</i>	customer
<i>Streptococcus pneumoniae</i>	customer
<i>Trametes spec.</i>	customer
<i>Vibrio fischerii</i>	MN

### Various applications

DNA was successfully isolated from the above mentioned microorganisms by the NucleoSpin® Microbial DNA kit.

www.mn-net.com

**MACHEREY-NAGEL**



MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

Germany and international:  
Tel.: +49 24 21 969-0  
Fax: +49 24 21 969-199  
E-mail: info@mn-net.com

Switzerland:  
MACHEREY-NAGEL AG  
Tel.: +41 62 388 55 00  
Fax: +41 62 388 55 05  
E-mail: sales-ch@mn-net.com

France:  
MACHEREY-NAGEL EURL  
Tel.: +33 388 68 22 68  
Fax: +33 388 51 76 88  
E-mail: sales-fr@mn-net.com

USA:  
MACHEREY-NAGEL Inc.  
Tel.: +1 484 821 0984  
Fax: +1 484 821 1272  
E-mail: sales-us@mn-net.com

